3.7.1 Camera Based Beam Near-Field Propagation Analyzer: M² 3.7.1.1 BeamSquared[®]

- ISO compliant
- · Automatically measure your beam quality in under 1 minutes
- Tune your laser for best operation
- Specifically developed for continuous usage
- Unequaled accuracy using patented Ultracal[™] Calibration
- Long optical train & automatic attenuation adjustment
- Flexible mounting configurations, install horizontal or vertically
- Pulsed and CW for most beam diameters and powers
- Compact and portable
- Detectors from 266nm to 10.6µm

The BeamSquared[®] system is a compact and fully automated tool for measuring the propagation characteristics of CW and pulsed laser systems from the UV to NIR to Telecom wavelengths. Users can also measure wavelengths above 1.8 microns, including CO₂ and terahertz in manual mode (a bench set-up; without the automated optical train) with a Pyrocam[™] IV or IIIHR. Our longer optical train and patented Ultracal[™] Calibration makes BeamSquared the most accurate product on the market and is ISO 11146 compliant. Its operational robustness and reliability ensures continuous use applications in industry, science, research and development.


Automatic M² - at Production Speeds


The BeamSquared optical train uses a fixed position lens with movable mirrors and camera. The mirrors that direct the focused beam into the camera are moved to precise locations, translating the beam through the near field, the waist, and the far field regions. All these measurements and translations, as well as incremental beam attenuation, are automatically controlled by the BeamSquared software. Design improvements in the BeamSquared system have decreased the measurement reporting time by 2-3 times, making it possible to report M² in under a minute.

Manual M²

Manual mode is available for wavelengths greater than NIR, particularly Terahertz and above, and for beams that are too large or too small for the BeamSquared optical system. Users are required to provide a manual translation/ attenuation apparatus.



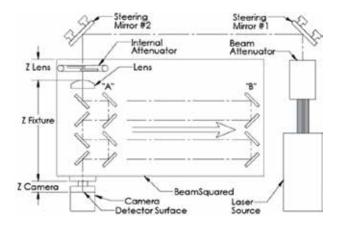

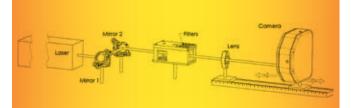

Software Features

Features	
Measurements	
	M ² x, M ² y, Kx, Ky, BPPx, BPPy
	Width at waist Wx, Wy
	Divergence angle Qx, Qy
	Waist location Zx, Zy
	Rayleigh X, Y
	Astigmatism
	Asymmetry ratio
	Statistical results are available on all measurements
Supports both automated a	nd manual runs
New Hardware	
	Allows standalone computer control of BeamSquared motion system
	Camera Options include: SP920, SP204S, Xeva, Pyrocam™ III HR or IV
	RF Lens Reader
	Lens must be present for operation
	Lens configuration data stored with lens (Focal length, calibration wavelength, material, etc.)
	Shutter only open what in live mode
	Table and attenuator calibration at startup (homing before each run)
Supports hardware Trigger	
Faster run times than M ² -20	05
New Interface	
	Selectable theme colors
	Splash screen with progress bar
2D display	Splash screen with progress bar
2D display	Selectable Color Palette
	Manual Cursor when not running (Cursor at centroid otherwise)
Caustic Display	Manual Cursor when not running (Cursor at centroid otherwise)
Caustic Display	Selecting individual frames
	Selecting individual frames
	Auto Aperture
Dura lafa Diantau	Exclude points from run
Run Info Display	
	Displays Caution Notice when beams are non-conforming: (too dark, too bright, misaligned, too large or too small)
	Option to ignore misaligned beams
	th, Laser to box distance, Laser to lens and focal length in manual mode)
Calculations	
	Frame Results (Total, Min, Peak, % in Aperture, Avg Pwr Density, Beam Width, Centroid, Peak, Cross Sectional Area)
	Laser Results (Waist Width, Divergence, Waist Location Rayleigh Length, M ² , K, BPP, Astigmatism, Asymmetry)
	After Lens Results (Waist Width, Divergence, Waist Location Rayleigh Length, Astigmatism, Asymmetry)
	Effective Focal Length of lens
	Fitted/Measured Divergence
	Supported Beam Width calculations
	• D4 Sigma
	Knife Edge 10/90 and Programmable
	EPSA - Encircled Power Smallest Aperture (power in a bucket)
Multiple Runs	
	Result statistics
	Progress Indicator
Single Page Report	
	Setup information
	Results
	Statistics
	Caustic chart
Logging/Export data	



Accuracy by Design

Spiricon products are known for accuracy. Using our patented Ultracal calibration method, auto aperturing to exclude noise beyond the wings of the laser beam, and long optical path, assures the user of the most accurate measurements in the industry.

Designed by Our Customers

Guided by customer input from our widely deployed previous generation M²-200s system, Spiricon redesigned the BeamSquared® to meet the challenging demands of the laser industry. The new BeamSquared system has significantly higher durability and operational robustness for continuous use in a three shifts a day, seven days a week environment. The rigid baseplate and internal optics greatly simplifies and reduces the time for initial set-up and alignment. The lens configuration data is now stored using an RF ID chip embedded in the lens holder which is uploaded automatically by the BeamSquared system when the lens cartridge is inserted in the system, eliminating the need for our customers to keep track of configuration file. Both novice and seasoned users will appreciate these new features along with the time-tested excellence that Spiricon has provided over the years.

Measurements


BeamSquared measures propagation characteristics in both the X and Y axes and displays the following parameters:

- Waist diameters
- Full angle Divergences
- Waist locations
- Rayleigh lengths
- M² or K and BPP
- factors
- Astigmatism
- Asymmetry

To optimize bench space, BeamSquared can be mounted either horizontally or vertically. Laser beam input port is the same dimension with either mounting method, X = Y, and the same as the M²-200s that it is replacing.



Main Screen Functions

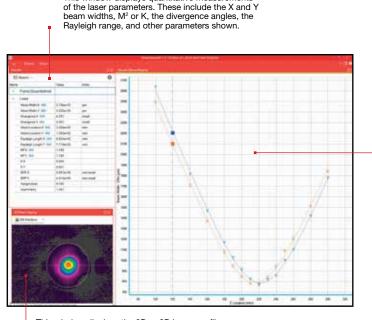

This window displays quantitative measurements

This window displays the 2D or 3D beam profile of the currently measured point in the beam propagation curve. This image enables visual intuitive verification of the beam profile behavior through focus. After each run the user can click any individual measured point and observe the beam profile. Outlying or anomalous points can be automatically or manually excluded from the curve fit calculations for more accurate results. This window presents measurements of beam width vs. position for a given run. After measuring a few points, the software extrapolates a curve fit. The Xs and Ys represent individual measurement points. The solid lines present the best fit hyperbola of the beam propagation equation to the measured points. The M² and other laser parameters are computed from the best fit hyperbola since it provides a smoothing of the data points.

Specifications

Model	BSQ-SP204S	BSQ-XC130	BSQ-A	BSQ-PY-M	
Beam Profiling					
Camera model	SP204S	XC130	BeamSquared [®] software,	Pyrocam [™] IIIHR or Pyrocam [™] IV	
Sensor type	Silicon CMOS	InGaAs CCD	software license, and optical	Software only, camera and optical	
Vavelengths	266 – 1100nm ⁽¹⁾	900 – 1700nm ⁽²⁾	train, no camera included	train not included.	
Active area	6.7mm x 5.6mm	9.6mm x 7.6mm	,	See individual camera data sheets	
Elements	2472 x 2064	320 x 256			
Effective pixel	2.74µm x 2.74µm	30µm x 30µm			
·	67 dB	68dB			
Dynamic range					
rame rate	37 fps	100 fps			
nterface	USB 3.0 ±5% typical, ±10% waist location and Rayleigh length typical				
Accuracy					
leasurement cycle time	<1 minute typical, depending				
Camera attachment		I C-mount, 90° camera on axis rotation			
ranslation system	Step-motor driven ball screw				
ranslation stage resolution	0.05mm				
tandard optics					
Lenses included (3)	266-440nm UV 500mm FL 430-700nm VIS 500mm FL 650-1000nm NIR 400mm FL 1000-1700nm Extended NIR 600mm FL	1000-1700nm Extended NIR 400mm FL 650-1000nm NIR 400 FL	266-440nm UV 500mm FL 430-700nm VIS 500mm FL 430-700nm VIS 400mm FL 650-1000nm NIR 400mm FL 1000-1700nm Extended NIR 400mm FL	N/A	
Optional lenses ^{(4) (5)} 'Not part of the standard BSQ kit	266-440nm UV 750mm FL 266-440nm UV 1000mm FL 430-700nm VIS 750mm FL 430-700nm VIS 1000mm FL 650-1000nm NIR 750mm FL 1000-1550nm Extended NIR 750mm FL	650-1000nm NIR 750mm FL 1000-1550nm Extended NIR 750mm FL	266-440nm UV 750mm FL 266-440nm UV 1000mm FL 430-700nm VIS 750mm FL 430-700nm VIS 1000mm FL 650-1000nm NIR 750mm FL 1000-1550nm Extended NIR 750mm FL	N/A	
Attenuation range					
(6)	Nominally from ND 1.0 to ND	4.8. Actual values vary with wa	velength.	N/A	
Damage limits ⁽⁶⁾					
	0.15 mW/cm ² CW mode 1.0 μ J/cm ² pulse mode Both of the above for an M ² =1 @ 1064nm	100 mW/cm ²	Depends on type of the camera	See camera data sheets	
Optical limits					
Vavelengths (7)	266 - 1100nm	900 - 1700nm	Depends on type of the camera	1.06 - 3000µm	
Beam size	BeamSquared Auto Mode 1m Varies with wavelength, waist			Pyrocam IIIHR 0.8mm – 10mm ma Pyrocam IV 0.8mm – 20mm max Depends on customer mechanics and lens	
Vinimum beam width	27µm	300µm	N/A	800µm	
Software					
BeamSquared Software	Fast scan method (1 minute) f	or automatic (ISO) and manual	M ² measurement, Windows 11 (6	64)	
General		, , , , , , , , , , , , , , , , , , , ,	,		
Storage temperature	-30° C to 65° C			N/A	
Storage humidity	95% maximum (non-condens	ing)		N/A	
Operating temperature	10° C to 40° C			N/A	
Operating humidity	95% maximum (non-condens	ing)		N/A	
Power requirements ⁽⁸⁾	55 % maximum (non-condens			IN/75	
	90 – 264 V AC			N/A	
nput voltage					
AC Line current	1.6 A			N/A	
ine frequency	47Hz to 63Hz			N/A	
Veight	26 lbs. w/o camera			N/A	
Dimensions	217.2mm X 459.5mm X 156.3	mm		N/A	
Compliance	CE, UKCA, China RoHS				
Ordering information					
Part Number	SP90630	SP90444 ⁽⁹⁾	SP90445	SP90410	

Notes:

 Number
 SP90630
 SP90444
 (9)
 SP9045
 SP90410

 (1) If device is used below 300 nm, measurement accuracy may degrade and long-term intensive irradiation at UV wavelengths may cause permanent damage to the imager.
 (2) For wavelengths between 1300-1400nm inner reflections have been observed that may impact beam measurement
 (3) Different lenses are required for different wavelength regions, spot sizes and divergences, Additional lenses must be ordered separately.
 (4) For selection of optimal BeamSquared lens, use Beam Profiler Finder.

 (5) These lenses have been calibrated for measurement of lasers with Rayleigh length up to 20 meters, for longer Rayleigh lengths, please use Beam Profiler Finder.
 (6) Damage threshold listed is based on products with similar design. Silicon cameras can be damaged by power in excess of 0.15 mW/cm² or energy in excess of 1 µJ/cm². BeamSquared employs a focusing optic. While it may be that the laser input power or energy measures well below this damage threshold, it can easily exceed these levels when focused onto the camera sensor. Use caution and error on the side of safety. Cameras can be costly to repair or replace.
 (7) For UV lasers 245-440nm, it is recommended to use the UV reflective ND2 filter P/N SP90568

 (8) For the optical train only. The PC computer supplies the power for the system components, such as the USB3 camera.
 (9) P/N for USA only, for any other country please consult Ophir representative





Accessories Ordering Information

Item	Description	P/N
BSQ-SP204S-A	An SP204S camera licensed for BeamSquared [®] . Sold as an accessory for those also purchasing a BSQ-XC130	SP90631
BSQ-Lens Kit UV-XNIR	Lens kit that includes 4 BeamSquared lenses: 500mm UV, 500mm VIS, 400mm NIR, 600mm XNIR	SP98009
BSQ-Lens Kit 266-1550	Lens kit that includes 5 BeamSquared lenses: 500mm UV, 500mm VIS, 400mm VIS, 400mm NIR, 400mm XNIR	SP90449
BSQ-Lens Kit 650-1700	Lens kit that includes 2 BeamSquared lenses: 400mm NIR, and 400mm XNIR.	SP90450
BSQ-Lens UV 500mm	Single BeamSquared lens, 500mm focal length, A/R coated for 266-440nm	SP90451
BSQ-Lens VIS 500mm	Single BeamSquared lens, 500mm focal length, A/R coated for 430-700nm	SP90452
BSQ-Lens VIS 400mm	Single BeamSquared lens, 400mm focal length, A/R coated for 430-700nm	SP90453
BSQ-Lens NIR 400mm	Single BeamSquared lens, 400mm focal length, A/R coated for 650-1000nm	SP90454
BSQ-Lens XNIR 400mm	Single BeamSquared lens, 400mm focal length, A/R coated for 1000-1550nm	SP90455
BSQ-Lens XNIR 600mm	Single BeamSquared lens, 600mm focal length, A/R coated for 1000-1550nm	SP90485
BSQ-Lens UV 750mm	Single BeamSquared lens, 750mm focal length, A/R coated for 266-440nm	SP90554
BSQ-Lens VIS 750mm	Single BeamSquared lens, 750mm focal length, A/R coated for 430-700nm	SP90555
BSQ-Lens NIR 750mm	Single BeamSquared lens, 750mm focal length, A/R coated for 650-1000nm	SP90556
BSQ-Lens XNIR 750mm	Single BeamSquared lens, 750mm focal length, A/R coated for 1000-1550nm	SP90557
BSQ-Lens UV 1000mm	Single BeamSquared lens, 1000mm focal length, A/R coated for 266-440nm	SP90558
BSQ-Lens VIS 1000mm	Single BeamSquared lens, 1000mm focal length, A/R coated for 430-700nm	SP90559
BSQ SP300 or SP920 to SP204S Upgrade	Camera upgrade	SP90632
BGS license for BSQ-SP204S	Includes BeamGage Standard software license in addition to BeamSquared software license	SP90633
BGP license for BSQ-SP204S	Includes BeamGage Professional software license in addition to BeamSquared software license	SP90634
BGP license for BSQ-XC130	Includes BeamGage Professional software license in addition to BeamSquared software license	SP90508
BSQ-XC130-KEY	Includes BeamSquard software license for XC-130 camera	SP90503
BSQ - UV Reflective Filter	BeamSquared reflective ND2 filter, UV Grade Fused Silica, Inconel coating for 245-440nm	SP90568

